INSTRUCTION MANUAL

STEREO ZOOM MICROSCOPE XTS2021

Thank you for purchasing the Stereo Zoom Microscope. Please read this instruction manual thoroughly before use to insure proper operation and long service.


NOTES

- This is a precision instrument – Handle With Care. Use caution to avoid shock or excessive forces on instrument when in use or transport.
- Avoid direct sunlight or areas with high temperatures or high levels of dust. Keep in location where temperate and humidity are stable.
- Keep lens surface clean from dirt and fingerprints.
- Clean the lens gently with lens cloth or soft cotton cloth.
- Do not turn the left and right Zoom Controls in opposite directions; this will damage the instrument.
- When not in use store with protective cover, in location that has low moisture in order to prevent internal mold growth.
- Adjust the Light Power Dial to adjust the brightness of the epi-illuminator for reflected light. When first turned on, rotate the dial to check that the power to the light is on.
- Lighting unit not certified for use outside of Japan.

SPECIFICATIONS

Rated Voltage	AC100V 50/60Hz
Magnification	6.5~45× (for 10× eyepiece)
	13~90× (for 20× eyepiece)
Objective Lens	0.65~4.5× (Zoom Ratio 1:7)
Eyepiece Lens	WF10× (FOV = 20mm) • WF20× (FOV = 10mm)
Working Distance	95mm
Eyepiece Angle	45°
Focus Adjustment	Movable Body
Stand	Rack and Pinion Type
Stage Plate	φ95mm×2pcs ①Ground Glass ②black/white (plastic)
Eye Width Adj. Range	55~75mm
Diopter Adj. Range	±5diopter
Illumination	Epi-illuminator: 6V • 15W Halogen bulb
	Transmitted: 5W ring light
Base	L245×W190×H60mm
Weight	approx.5.7kg

PART IDENTIFICATION & SETUP

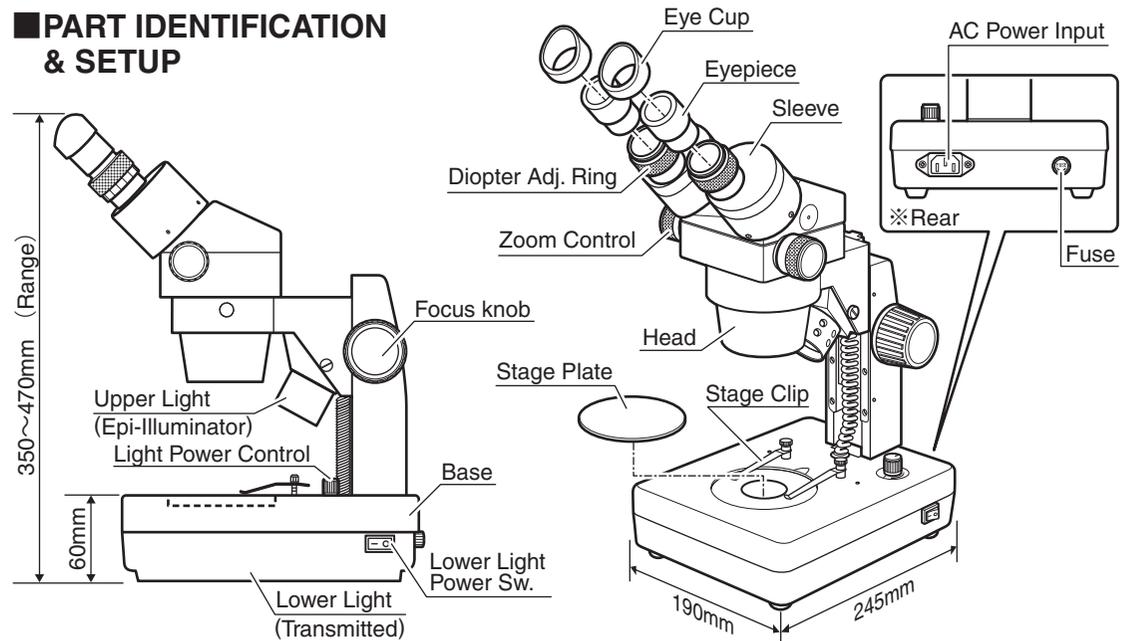OPERATION

SAMPLE PLACEMENT

- ① Place the sample in the center of the Stage Plate
- ② If required, use Stage Clips to hold sample in place.

EYE CUP USE

- For use without eyeglasses : Hold Eyepiece and rotate the Diopter Adjustment Ring to adjust diopter setting. Align the Eye cups to point toward user and look through the Eyepiece with eyes brought close.
- For use with eyeglasses : Remove Eye cups to use.

EYE WIDTH ADJ.

Hold both sleeves in each hand and look through the Eyepieces. Adjust the Sleeves to align the left and right field of view seen through the Eyepieces.

FOCUS & DIOPTR ADJUSTMENT

- ① Rotate the Zoom Control for highest magnification.
- ② Look through the right eyepiece with the right eye and adjust the focus using the Focus Knob.
- ③ Rotate the Zoom Knob for lowest magnification.
- ④ Using the Right Diopter adjustment, adjust the focus if required.
- ⑤ Now rotate the Zoom Control back to highest magnification and check the focus. Repeat steps ② ~ ④ to achieve the best focus accuracy.
- ⑥ Rotate the Zoom Control for highest magnification. Look through the left Eyepiece with the left eye, and adjust focus using the Left Diopter Adjustment Ring.

FUSE REPLACEMENT

In case of fuse failure, it can be replaced with a commercially available 250V 2A fuse.

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