

LED LIGHT SCALE LOUPE LSL-26/LSL-35

Thank you for purchasing the Led Light Scale Loupe.
Please read this manual thoroughly before use to insure proper operation and long service.

FEATURES

- Magnifying glass with scale and compact LED light

ACCESSORIES

- Scale Reticle, 2× (1× installed)
- Lens Cloth, 1×
- Eyepiece Cap, 1×

SPECIFICATIONS

Model No.	Magnification	Field of View	Effective Diameter	Reticle Diameter	LED Light Continuous On Time	Power Supply
LSL-26	10×	24mm	16.5mm	26mm	approx. 10hr. (for new battery)	LR41×3 (sold separately)
LSL-35	10×	33mm	20mm	35mm		

PRODUCT NOTES


CAUTION

- Never look at sun or strong light through the Loupe, there is risk of eye damage.
- Do not expose skin to sunlight or strong light through the Loupe, there is risk of burns.
- Do not leave Loupe in direct or indirect sunlight, there is risk of fire.

NOTICE


- To prevent damage, do not drop or mishandle.
- Protect from high temperatures and humidity to prevent degradation and damage.
- Clean lens with lens cloth and use gentle motion to prevent dirt from scratching surface.
- Remove protective film before use.
- When Scale Reticle is not in use, store in protective case.
- Use only as directed.

PARTS IDENTIFICATION

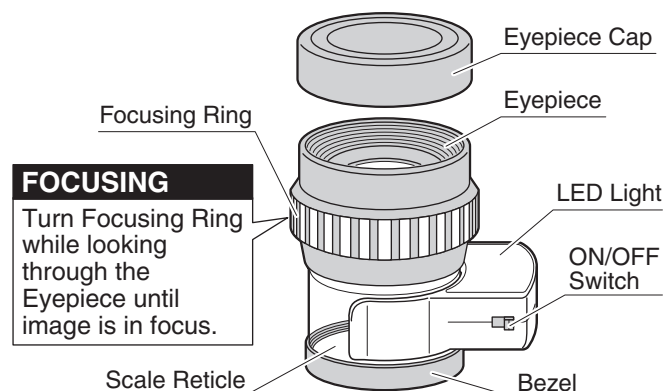

LED LIGHT


Mounting

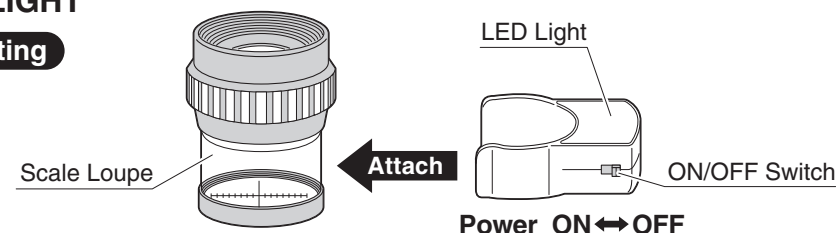

Battery Replacement

- ① Slide Battery Cover in direction of arrow.
- ② Insert 3 LR41 batteries, using care to put ⊕, ⊖ sides in proper direction.
- ③ Replace Battery Cover.

NOTICE Use only LR41 type battery.

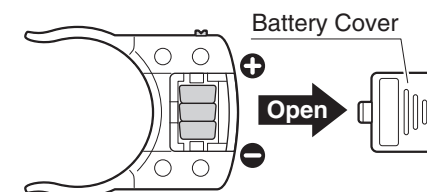


SCALE


Exchanging Scales

LSL-26

- ① Remove the Reticle Bezel and Reticle from Loupe.
- ② Replace the with the Reticle containing the desired Scale.

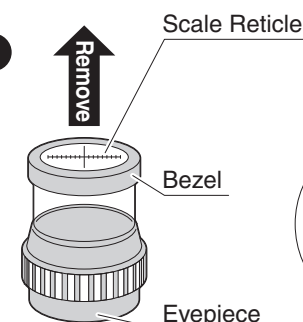

※ LSL-26 Reticle and Bezel are a single unit, please replace as pair without removing Reticle from Bezel.

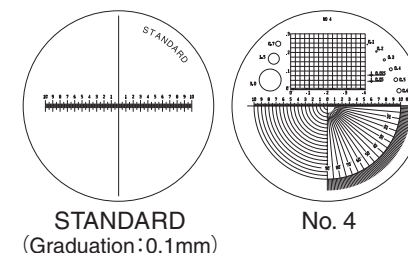

LSL-35

- ① Rotate the Bezel to remove.
- ② Place the desired Reticle on the Loupe.
- ③ Replace the Bezel and rotate to hold Reticle in place.

